


Datasheet
LCA-S-400K-BSI
**Low Noise 400 kHz Photoreceiver
with Black Silicon Photodiode**


Features	<ul style="list-style-type: none"> • Large area black silicon photodiode, 2 × 2 mm active diameter • Quantum efficiency approx. 90 % from 240 nm to 1000 nm • Bandwidth DC – 400 kHz • Amplifier transimpedance gain 1.0×10^7 V/A • Max. conversion gain 7.3×10^8 V/W @ 1010 nm • Spectral range 200 – 1100 nm • Free-space input 1.035"-40 threaded • UNC 8-32 and M4 tapped holes for mounting on standard posts with metric and imperial thread
----------	--

Applications	<ul style="list-style-type: none"> • Spectroscopy • General purpose opto-electronic measurements • Optical front-end for oscilloscopes, A/D converters and lock-in amplifiers
--------------	--



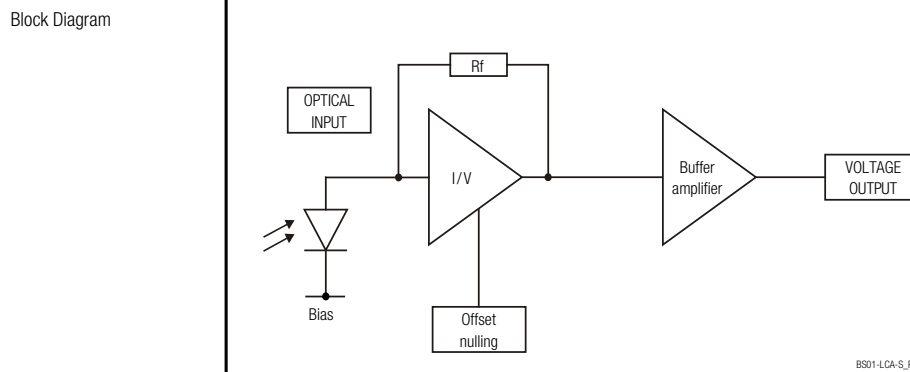





Intended Use	<p>The LCA-S-400K-BSI photoreceiver consists of a black silicon photodiode and a subsequent low-noise fixed gain transimpedance amplifier. It is designed for fast conversion of small optical signals into equivalent output voltages. Operation is mostly self-explanatory. If in doubt, consult this document or contact support@femto.de.</p> <p>For safe operation, please refer to the damage thresholds specified in the "Absolute Maximum Ratings", "Temperature Range" and "Power Supply" sections of this document.</p> <p>The operating environment must be free of smoke, dust, grease, oil, condensing moisture, and other contaminants that could affect the operation or performance.</p>
--------------	--

SOPHISTICATED TOOLS FOR SIGNAL RECOVERY
F E M T O

LCA-S-400K-BSI_R2/TH/19MAR2026

Datasheet
LCA-S-400K-BSI
**Low Noise 400 kHz Photoreceiver
with Black Silicon Photodiode**

Available Version	<p>LCA-S-400K-BSI-FST</p>  <p>1.035"-40 threaded flange with internally threaded coupler ring (outer diameter 30 mm) for free space applications, compatible with many optical standard accessories</p>
Related Models	<p>LCA-S-400K-SI-FST Si-PIN, Ø 3 mm, 320 - 1060 nm free space input, 1.035"-40 threaded flange</p> <p>LCA-S-400K-IN-FST InGaAs-PIN, Ø 0.5 mm, 900 - 1700 nm free space input, 1.035"-40 threaded flange</p>
Available Accessories	<p>PRA-FC  Fiber-adaptor with external PRA-FCA  1.035"-40 thread PRA-FSMA </p> <p>PRA-PAP  Alternative mounting option: post adapter plate, easy to mount on FEMTO photoreceiver series OE, FWPR, PWPR, HCA-S and LCA-S</p> <p>PS-15-25-L  Power Supply input: 100 – 240 VAC output: ±15 VDC</p>
Specifications	<p>Test conditions $V_s = \pm 15 \text{ V}$, $T_A = 25 \text{ }^\circ\text{C}$, output load impedance 1 MΩ, warm-up 20 minutes (min. 10 minutes recommended)</p> <p>Gain Transimpedance gain $1.0 \times 10^7 \text{ V/A}$ (@ output load $\geq 100 \text{ k}\Omega$) Gain accuracy $\pm 1 \%$ (electrical) Conversion gain $7.3 \times 10^6 \text{ V/W typ.}$ (@ 1010 nm, output load $\geq 100 \text{ k}\Omega$)</p> <p>Frequency Response Lower cut-off frequency DC Upper cut-off frequency (-3 dB) 400 kHz Gain flatness $\pm 0.5 \text{ dB}$</p> <p>Time Response Rise/fall time (10 % – 90 %) 900 ns</p> <p>Input Noise equivalent power (NEP) 110 fW/$\sqrt{\text{Hz}}$ (@ 1010 nm, 10 kHz) Optical saturation power 1.37 μW (for linear amplification, @ 1010 nm) Input offset compensation range $\pm 300 \text{ nA}$, adjustable by offset potentiometer</p>


SOPHISTICATED TOOLS FOR SIGNAL RECOVERY
F E M T O

LCA-S-400K-BSI_R2/TH/19MAR2026

Datasheet
LCA-S-400K-BSI
**Low Noise 400 kHz Photoreceiver
with Black Silicon Photodiode**

Specifications (continued)		
Detector	Detector Active area Spectral range Max. sensitivity	Black silicon photodiode 2 × 2 mm 200 – 1100 nm 0.73 A/W typ. (@ 1010 nm)
Output	Output voltage range Output impedance Max. output current Output noise	–3 V ... +10 V (@ ≥ 100 kΩ output load) 50 Ω (terminate with ≥ 100 kΩ load) 30 mA (short-circuit proof) 2.0 mV RMS (18 mV peak-peak) typ. (@ ≥ 100 kΩ load, no signal on detector, measurement bandwidth 20 MHz)
Input Flange	Material	1.4305 stainless steel, nickel-plated
Coupler Ring	Material	1.4305 stainless steel, glass bead blasted
Power Supply	Supply voltage Supply current	±15 V (±14.5 V ... ±16.5 V) ±40 mA (depends on operating conditions, recommended power supply capability min. ±150 mA)
Case	Weight Material	212 g (0.47 lbs) incl. coupler ring AlMg4.5Mn, nickel-plated
Temperature Range	Storage temperature Operating temperature	–30 °C ... +85 °C 0 °C ... +60 °C

Absolute Maximum Ratings	Optical input power (CW) Power supply voltage	10 mW ±20 V
--------------------------	--	----------------

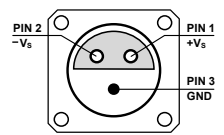Connectors	Input Output Power supply	1.035"-40 threaded flange for free space applications and for use with various types of optical standard accessories BNC jack (female) LEMO® series 1S, 3-pin fixed socket (mating plug type: FFA.1S.303.CLAC52)
		 <p>PIN 1 +Vs PIN 2 -Vs PIN 3 GND</p> <p>Pin 1: +15 V Pin 2: –15 V Pin 3: GND</p>

Scope of Delivery	LCA-S-400K-BSI, internally threaded coupler ring, LEMO® 3-pin connector, datasheet, transport package
-------------------	---


Ordering Information	LCA-S-400K-BSI-FST	Black silicon photodiode, 2 × 2 mm, 200 – 1100 nm free space input, 1.035"-40 threaded flange
----------------------	--------------------	--

SOPHISTICATED TOOLS FOR SIGNAL RECOVERY
F E M T O


LCA-S-400K-BSI_R2/TH/19MAR2026

Datasheet
LCA-S-400K-BSI
**Low Noise 400 kHz Photoreceiver
with Black Silicon Photodiode**

Spectral Responsivity

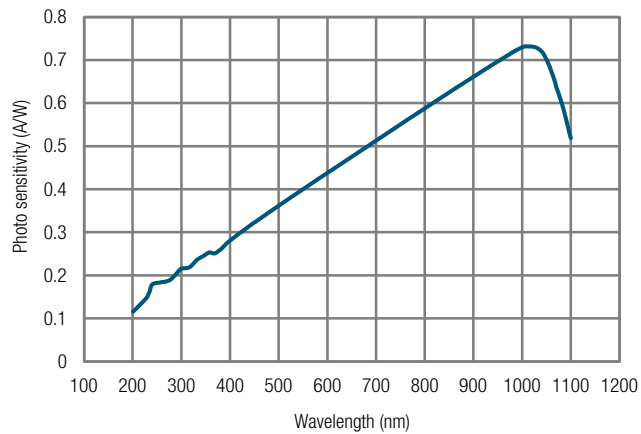

Spectral Quantum Efficiency

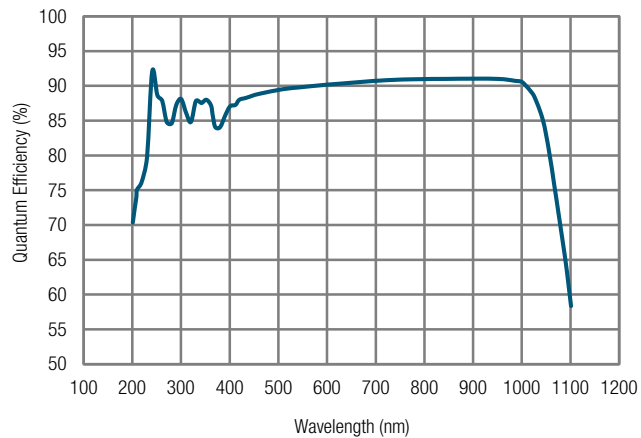

SOPHISTICATED TOOLS FOR SIGNAL RECOVERY


F E M T O

LCA-S-400K-BSI_R2/TH/19MAR2026

Datasheet
LCA-S-400K-BSI
**Low Noise 400 kHz Photoreceiver
with Black Silicon Photodiode**

Dimensions

LCA-S-400K-BSI-FST (1.035"-40 threaded free space input)

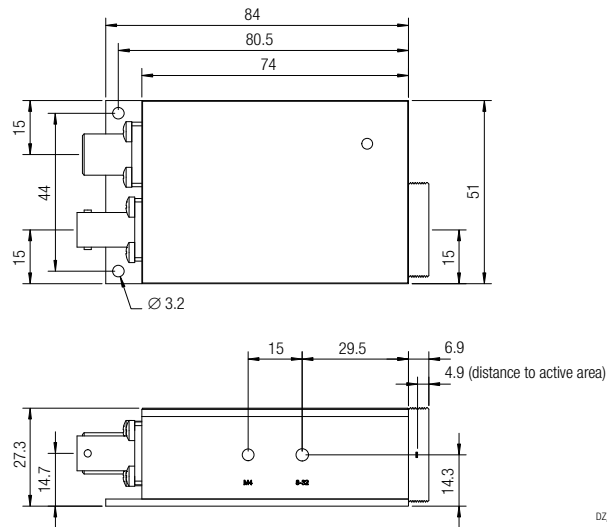

DZ_LCA-S-400K_R1

all dimensions in mm unless otherwise noted

Specifications are subject to change without notice. Information provided herein is believed to be accurate and reliable. However, no responsibility is assumed by FEMTO Messtechnik GmbH for its use, nor for any infringement of patents or other rights of third parties which may result from its use. No license is granted by implication or otherwise under any patent or patent rights of FEMTO Messtechnik GmbH. Product names mentioned may also be trademarks used here for identification purposes only.

© by FEMTO Messtechnik GmbH · Printed in Germany

SOPHISTICATED TOOLS FOR SIGNAL RECOVERY

F E M T O

LCA-S-400K-BSI_R2/TH/19MAR2026